



ELECTRONIC BALANCES AND SCALES

SUBSTANCE – SCRUBBING MILK – LEMON

Below information on determination of humidity content for above specified substance is a supplement of traditional drying method. RADWAG technical support department hopes, that this datasheet will be helpful for the tests performed with application of RADWAG moisture analyzers.

DRYING PARAMETERS – TRADITIONAL METHOD

Substance description:	Scrubbing milk – lemon
Reference /traditional/ method parameters:	105 °C / 4 hours.
Dry masscontent determined with traditional method:	53,89 %
Standard deviation:	0,025 %



DRYING PARAMETERS - MOISTURE ANALZYER SERIES MAC 50

Equipment necessary for drying process	Disposable aluminum pans
Sample preparation	Place a thin layer over all pan surface
Sample size	~ 2,2 g
Drying profile	Standard
Drying temeprature	125 °C
Auto switch-off criterion	Automatic 3 [Auto 1 - 1mg/10sek; Auto 2 - 1mg/25sek; Auto 3 - 1mg/60sek; Auto 4 - 1mg/90sek; Auto 5 - 1mg/120sek]
Average measurement time	~ 12 minutes
Determined dry mass content /average value/	53,81 %
Standard deviation	0,097 %

© 2008 RADWAG. All Rights Reserved.

All copyrights to the materials contained in the publication belong to RADWAG and may not be copied, processed or used for commercial purposes without their knowledge or consent RADWAG.